

# Instructions for making reagents

## **Complete RPMI (cRPMI)**

Supplement sterile RPMI-1640 medium with 10% sterile heat-inactivated FBS and 1% sterile antibiotic/antimycotic. Store at 4°C.

## **Stock 25% Human Serum Albumin (HSA)**

Prepare 25% has in RPMI; e.g. dissolve 25gm human albumin fraction V in 100ml RPMI (**not** complete RPMI). Allow 24 hours to fully dissolve. Store at 4°C.

## **12.5% Human Serum Albumin (HSA)**

Combine 10ml of stock 25% HAS and 10ml of sterile RPMI (**not** complete RPMI)

## **2x Freezing Medium**

Combine 10ml of stock 25% human serum albumin (HSA) and 10ml of sterile RPMI (**not** complete RPMI). Add 5ml DMSO. Store at 4°C.